

WEB INSERT

130113



**ARBOR
ASSAYS**

DetectX[®]
Ceruloplasmin
Colorimetric Activity Kit

2 Plate Kit

Catalog Number K035-H1

SPECIES INDEPENDENT

Sample Types Validated:

Serum and Urine samples

Protected by US Patent 8,227,206

Please read this insert completely prior to using the product.

FOR RESEARCH USE ONLY. NOT FOR USE IN DIAGNOSTIC PROCEDURES.

www.ArborAssays.com

WEB INSERT
TABLE OF CONTENTS

130113

Background	3
Assay Principle	4
Related Products	4
Supplied Components	5
Storage Instructions	5
Other Materials Required	6
Precautions	6
Sample Types	7
Sample Preparation	7
Standard Preparation	8
Assay Protocol	9
Calculation of Results	10
Typical Data	10-11
Validation Data Sensitivity, Linearity, etc.	11-13
Sample Values	14
Warranty & Contact Information	15
Plate Layout Sheet	16



Ceruloplasmin (Cp) is a multicopper oxidase enzyme involved in the safe handling of oxygen in some metabolic pathways of vertebrates. Discovered in 1948, a blue protein from the α_2 -globulin fraction of human serum possessing oxidase activity towards aromatic diamines and catechol was purified by Holmberg and Laurell¹. It was denoted ceruloplasmin, literally meaning 'a blue substance from plasma'. Specialized copper sites have been recruited during evolution to provide long-range electron transfer reactivity and oxygen binding and activation in proteins destined to cope with oxygen reactivity in different organisms. Ceruloplasmin belongs to the family of multi-copper oxidases which are among the few enzymes able to bind molecular oxygen to perform its complete reduction to water^{2,3}. Ceruloplasmin contains 95% of the copper in serum⁴. Cp found in serum is expressed in the liver, but it is also expressed in the brain, lung, spleen and testis.

Aceruloplasminaemia is an autosomal recessive disorder of iron metabolism characterized by the complete absence of ceruloplasmin^{5,6}. The role of Cp in tissue iron overload and the subsequent clinical findings of diabetes, retinal degeneration and neurodegeneration has been associated with iron overload in aceruloplasminaemic patients⁷. Thus it is clearly indicated that ceruloplasmin plays an essential role in iron metabolism. Ceruloplasmin is also associated with reproduction. Copper-deficient female rats seem to be protected against mortality. This protection has been suggested to be provided by estrogens, since estrogens alter the subcellular distribution of copper in the liver, an increase in plasma copper levels and subsequent ceruloplasmin synthesis⁸.

1. Holmberg CG and Laurell CB. Investigation serum copper. II. Isolation of the copper containing protein, and a description of some its properties. 1948, Acta. Chem. Scand. 2:550-556.
2. Malmstrom BG. Enzymology of oxygen. Annu. Rev. Biochem. 1982, 51: 21-59.
3. Farver O. and Pecht I. Electron transfer reactions in multi-copper oxidases. In: Multi-Copper Oxidases, 1997, pp. 355-389, Messerschmidt A. (ed.), World Scientific Publication, Singapore.
4. Harris, ZL and Gitlin, JD, Genetic and molecular basis for copper toxicity, Am. J. Clin. Nutr. 1996, 63:836S-841S.
5. Yoshida K., . et al. A mutation in the ceruloplasmin gene is associated with systemic hemosiderosis in humans. Nat. Genet. 1995, 9:267-272.
6. Harris ZL., Takahashi Y., Miyajima H., Serizawa M., MacGillivray RTA. and Gitlin JD. Aceruloplasminemia: molecular characterization of a novel disorder of iron metabolism. Proc. Natl. Acad. Sci. USA, 1995, 92:2539-2543.
7. Takahashi Y., Miyajima H., Shirabe S., Nagataki S., Suenaga, A. and Gitlin JD. Characterization of a nonsense mutation in the ceruloplasmin gene resulting in diabetes and neurodegenerative disease. 1996, Hum. Mol. Genet. 5:81-84.
8. Fields, M., Lewis, C., Scholfield, D., Powel, AS., Rose, AJ., Reiser, S., and Smith, JC., Female rats are protected against the fructose-induced mortality of copper deficiency. Proc. Soc. Expl Biol. Med. 1986, 186:145-149.

WEB INSERT ASSAY PRINCIPLE

130113

The DetectX[®] Ceruloplasmin Activity Kit is designed to quantitatively measure ceruloplasmin activity in diluted serum and urine samples. Please read the complete kit insert before performing this assay. A human ceruloplasmin standard is provided to generate a standard curve for the assay and all samples should be read off of the standard curve. Samples are diluted in the provided Assay Buffer and added to the wells of a half area clear plate. The reconstituted Ceruloplasmin Substrate is added and the plate is incubated at 30°C for 60 minutes. The ceruloplasmin in the standards and samples reacts with the substrate to produce a colored product. The optical density is read at 560 nm. Increasing levels of ceruloplasmin in the samples causes an increase in the fuschia (pink-purple) product. The activity of the ceruloplasmin in the sample is calculated after making a suitable correction for any dilution, using software available with most plate readers. The results are expressed in terms of units of ceruloplasmin activity per mL.

RELATED PRODUCTS

DetectX[®] Kits

Hemoglobin Dual Range Detection Kit	Catalog Number K013-H1
17 β -Estradiol EIA Kits	Catalog Number K030-H1/H5
Estrone EIA Kits	Catalog Number K031-H1/H5
Progesterone EIA Kits	Catalog Number K025-H1/H5
PGFM (13,14-dihydro-15-keto-Prostaglandin F _{2α}) EIA Kits	Catalog Number K022-H1/H5
Nitric Oxide Colorimetric Detection Kit	Catalog Number K023-H1
Superoxide Dismutase (SOD) Activity Kit	Catalog Number K028-H1

WEB INSERT
SUPPLIED COMPONENTS

130113

Clear 96 well Half Area Plates	2 Plates	Catalog Number X018-2EA
Ceruloplasmin Standard 1,000 Units/mL of human ceruloplasmin in a special stabilizing solution.	20 μ L	Catalog Number C120-20UL
Assay Buffer Concentrate A 5X concentrate containing detergents and stabilizers.	28 mL	Catalog Number X111-28ML
Ceruloplasmin Colorimetric Substrate Ceruloplasmin substrate lyophilized from a special stabilizing solution.	2 Vials	Catalog Number C121-1EA
Plate Sealers	2 Each	Catalog Number X002-1EA

STORAGE INSTRUCTIONS

This kit should be stored at -20°C until the expiration date of the kit.

Once opened the kit can be stored at 4°C up to the expiration date on the kit label, **except for the Ceruloplasmin Standard and reconstituted Ceruloplasmin Substrate, which must be stored at -20°C.**

WEB INSERT
OTHER MATERIALS REQUIRED

130113

Repeater pipet with disposable tips capable of dispensing 25 μ L.

An incubator capable of maintaining 30°C

96 well plate reader capable of reading optical density at 560 nm.

Software for converting optical density readings from the plate reader and carrying out four parameter logistic curve (4PLC) fitting. Contact your plate reader manufacturer for details.

PRECAUTIONS

As with all such products, this kit should only be used by qualified personnel who have had laboratory safety instruction. The complete insert should be read and understood before attempting to use the product.

The Ceruloplasmin Standard supplied in this kit is purified from human blood. The source of the protein was tested and found to be negative for hepatitis and HIV. However please treat the standard as a potentially infectious sample.

WEB INSERT
SAMPLE TYPES AND PREPARATION

130113

Samples that need to be stored after collection should be stored at -70°C or lower, preferably after being frozen in liquid nitrogen. This assay has been validated for serum and urine samples. Samples containing visible particulate should be centrifuged prior to using.

Ceruloplasmins are ancient enzymes that should behave in a similar manner to the colorimetric substrate. It is believed that the assay will measure Cp activity from a wide range of sources. It is up to the end user to determine if their samples can be measured using this assay.

SAMPLE PREPARATION

Serum and Urine Samples

Serum and urine samples should be diluted at least 1:20 fold in the Assay Buffer supplied.

For reporting urinary Ceruloplasmin Activity values use our Urinary Creatinine Detection kits, K002-H1 and K002-H5, to normalize to urine volume.

WEB INSERT

STANDARD PREPARATION

130113


Allow the kit reagents to come to room temperature for 30-60 minutes. We recommend that all standards and samples be run in duplicate to allow the end user to accurately determine Ceruloplasmin activities. Ensure that all samples have reached room temperature and have been diluted as appropriate prior to running them in the kit.

Assay Buffer

Dilute Assay Buffer Concentrate 1:5 by adding one part of the concentrate to four parts of deionized water. Once diluted this is stable at 4°C for 3 months.

Standard Preparation

Standards are prepared by labeling seven tubes as #1 through #7. Add 995 μL of Assay Buffer to tube #1. Pipet 300 μL of Assay Buffer into tubes #2 to #7. Carefully add 5 μL of the Ceruloplasmin Stock from the vial to tube #1 and vortex completely. Take 600 μL of the Cp solution in tube #1 and add it to tube #2 and vortex completely. Repeat the serial dilutions for tubes #3 through #7. The ceruloplasmin activity in tubes 1 through 8 will be 5, 3.33, 2.22, 1.48, 0.988, 0.658, and 0.439 U/mL.

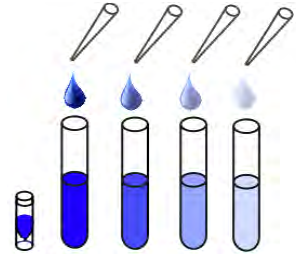

Use all Standards within 2 hours of preparation.

	Std 1	Std 2	Std 3	Std 4	Std 5	Std 6	Std 7
Assay Buffer Vol (μL)	995	300	300	300	300	300	300
Addition	Stock	Std 1	Std 2	Std 3	Std 4	Std 5	Std 6
Vol of Addition (μL)	5	600	600	600	600	600	600
Final Activity (U/mL)	5.0	3.33	2.22	1.48	0.988	0.658	0.439

Ceruloplasmin Substrate Preparation

Add 3 mL of water to the vial and mix thoroughly. This solution can be stored at 4°C for up to 2 weeks. The solution can also be stored at -20°C for up to the expiration date on the kit label.

WEB INSERT ASSAY PROTOCOL

130113

Use the plate layout sheet on the back page to aid in proper sample and standard identification. Set plate parameters for a 96-well Corning Costar 3695 plate. See: <http://www.ArborAssays.com/resources/lit.asp> for plate dimension data.

1. Pre-warm incubator to 30°C.
2. Pipet 100 μ L of diluted samples or appropriate standards into duplicate wells in the plate.
3. Pipet 100 μ L of Assay Buffer into duplicate wells as the Zero standard.
4. Add 25 μ L of the reconstituted Cp Substrate solution to each well using a repeater pipet.
5. Incubate at 30°C for 60 minutes.
6. Read the optical density generated from each well in a plate reader capable of reading at 560 nm.

WEB INSERT
CALCULATION OF RESULTS

130113

Average the duplicate OD readings for each standard and sample. Create a standard curve by reducing the data using the 4PLC fitting routine on the plate reader, after subtracting the mean OD for the zero standard. The sample activity obtained should be multiplied by the dilution factor to obtain neat sample values.

TYPICAL DATA

Sample	Mean Net OD	Ceruloplasmin Activity (U/mL)
Standard 1	0.922	5
Standard 2	0.663	3.33
Standard 3	0.428	2.22
Standard 4	0.290	1.48
Standard 5	0.152	0.988
Standard 6	0.108	0.658
Standard 7	0.049	0.439
Zero	0.000	0
Sample 1	0.337	1.831
Sample 2	0.209	1.169


**Always run your own standard curves for calculation of results.
Do not use these data.**

Ceruloplasmin Unit Definition

One Unit of Ceruloplasmin causes an increase in OD of 0.01 per minute at 37°C and pH 5.5 using N,N-dimethyl-p-phenylene diamine as substrate.



Typical Standard Curve

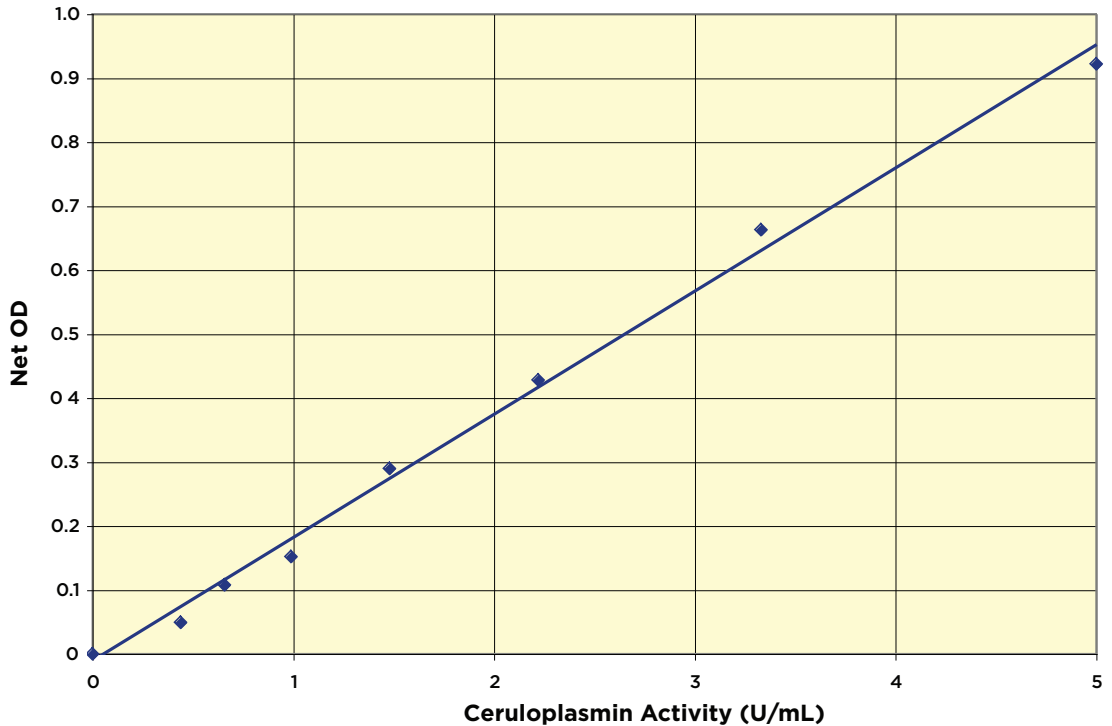

**Always run your own standard curves for calculation of results.
Do not use these data.**

VALIDATION DATA

Sensitivity

Sensitivity was calculated by comparing the ODs for twenty wells run for each of the zero and standard #7. The detection limit was determined at two (2) standard deviations from the zero along the standard curve.

Sensitivity was determined as 0.242 U/mL. This is equivalent to 24.2 mU/well.

Limit of Detection

The Limit of Detection for the assay was determined in a similar manner by comparing the ODs for twenty runs for each of the zero standard and a low concentration panda urine sample.


Limit of Detection was determined as 0.425 mU/mL. This is equivalent to 42.5 mU/well.

Linearity


Linearity was determined by taking two diluted panda urine or human serum samples, one with a high known ceruloplasmin activity and the other with a lower ceruloplasmin activity and mixing them in the ratios given below. The measured activities were compared to the expected values based on the ratios used.

High Sample	Low sample	Expected Activity (U/mL)		Observed Activity (U/mL)		% Recovery	
		Urine	Serum	Urine	Serum	Urine	Serum
80%	20%	2.10	1.40	2.22	1.27	105.5	90.5
60%	40%	1.72	1.54	1.77	1.71	102.6	111.0
40%	60%	1.35	1.68	1.57	1.67	116.5	99.6
20%	80%	0.97	1.82	0.97	1.99	99.5	109.5
Mean Recovery						106.0%	102.6%

Urine Linearity

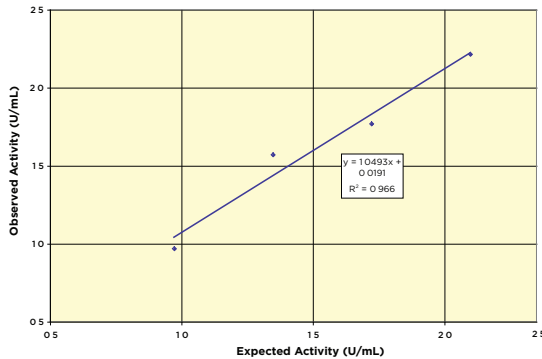

Serum Linearity

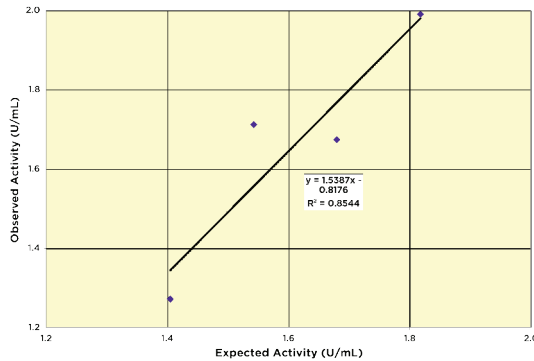

Intra Assay Precision

Two panda urine samples diluted in Assay Buffer were run in replicates of 20 in an assay. The mean and precision of the calculated activities were:

Sample	Ceruloplasmin Activity (U/mL)	%CV
1	1.85	5.0
2	1.17	7.1

Inter Assay Precision

Two panda urine samples diluted in Assay Buffer were run in duplicates in sixteen assays run over multiple days by three operators. The mean and precision of the calculated activities were:

Sample	Ceruloplasmin Activity (U/mL)	%CV
1	1.92	8.7
2	1.17	12.2

WEB INSERT SAMPLE VALUES

130113

Urine samples from a variety of mammals, some of them pregnant, were tested in the assay. Activity values, after adjustment for dilution, ranged from 11.26 U/mL to 73.7 U/mL, with an average of 29.6 U/mL. After adjusting for urinary creatinine the normalized activity values ranged from 11.3 to over 180 U/mg creatinine. Urinary Creatinine concentrations were determined using the Arbor Assays DetectX® kits K002-H1 and K002-H5.

This assay has not yet been tested on fecal extracts or plasma samples.

Seven random human serum were diluted with Assays Buffer and tested in the assay. Activity values ranged from 57.01 to 128.4 U/mL with an average value of 87.9 U/mL.

WEB INSERT
LIMITED WARRANTY

130113

Arbor Assays warrants that at the time of shipment this product is free from defects in materials and workmanship. This warranty is in lieu of any other warranty expressed or implied, including but not limited to, any implied warranty of merchantability or fitness for a particular purpose.

We must be notified of any breach of this warranty within 48 hours of receipt of the product. No claim shall be honored if we are not notified within this time period, or if the product has been stored in any way other than outlined in this publication. The sole and exclusive remedy of the customer for any liability based upon this warranty is limited to the replacement of the product, or refund of the invoice price of the goods.

CONTACT INFORMATION

For details concerning this kit or to order any of our products please contact us:

Arbor Assays

1514 Eisenhower Place
Ann Arbor, Michigan 48108 USA

Phone: 734-677-1774

Fax: 734-677-6860

Web: www.ArborAssays.com

E Mail Addresses:

Info@ArborAssays.com

Orders@ArborAssays.com

Technical@ArborAssays.com

Contracts@ArborAssays.com

DetectX®, ThioStar® and the Arbor Assays logo are all registered trademarks.

	A	B	C	D	E	F	G	H
1								
2								
3								
4								
5								
6								
7								
8								
9								
10								
11								
12								